

SEB7 extraction buffer (10X)

Catalog number: ACC 00505

Lot number	Item
_____	500 ml - makes 5 liter

Store SEB7 extraction buffer (10X) concentrate at room temperature.

Preparing buffer from concentrate

SEB7 extraction buffer is used for GMO composite seed testing.

Dilute 1 volume of 10X SEB7 extraction buffer concentrate with 9 volumes of distilled water before use.

Example: To prepare 2.5 liter of working SEB7 extraction buffer

Buffer concentrate	250 ml
Distilled water	2250 ml

Measure the distilled water into a container and add the buffer concentrate while stirring.

Storage

Agdia recommends preparing only as much 1X SEB7 extraction buffer as is needed for one day.