



## Intended Use:

This ImmunoStrip test is a rapid means of screening cucurbits and bacterial samples for *Acidovorax avenae* subsp. *citrulli* (Aac), the causal agent of bacterial fruit blotch. The Aac ImmunoStrip is intended for use with fruit, leaves, and seedlings. ImmunoStrip tests require no equipment or expertise to run. Results can be obtained in 30 minutes or less, making ImmunoStrips perfect for use in the field or greenhouse.

## Kit Storage:

Kit components should be stored refrigerated (2 - 8 °C) between uses, and ImmunoStrips should be tightly sealed in the desiccated container at all times.

Before use, allow all kit components to warm to room temperature (18 - 30 °C).

## Limitations:

Extraction and testing of decomposed, dried, or large amounts of tissue can cause invalid results.

### Contents of Kit:

- ImmunoStrips
- \*SEB4 sample extraction bags

### Not Included but Required:

- Scissors, knife, or razorblade
- Extraction tool
  - » Agdia tissue homogenizer ([ACC 00900](#)), marker, or pen
- Letter holder or another device to hold sample extraction bags upright
- \*\*1.5 mL microcentrifuge tubes
- \*\*Pipettes (100 µL & 1000 µL)

\*Not included if ordering STX only  
 \*\*Only for bacterial samples

## PERFORMING THE ASSAY (\*Special Attention Required)

### Prepare Sample

**1.** Samples should be taken from symptomatic tissue when possible. Agdia sample extract bags **contain 3 mL** of extraction buffer. Use the chart to the right to determine the sample size to achieve the optimal dilution for each sample type.

| Sample Type | Sample Size   |
|-------------|---------------|
| Seedling    | 0.15 g (1:20) |
| Leaf/Stem   | 0.15 g (1:20) |
| Early Fruit | 0.3 g (1:10)  |
| Late Fruit  | 0.6 g (1:5)   |

- **Seedling samples:** Symptoms include small water-soaked lesions that later become necrotic on the underside of cotyledons or leaves. (Figure 1)
- **Plant samples:** Symptoms include light brown to reddish brown lesions along the midrib of the leaf. (Figure 1)
- **Fruit samples:** For **early infection** symptoms include a small water-soaked area that expand with irregular margins. The sample should be wedge-shaped deep enough to get past rotten material. (Figure 2) For **late infection** symptoms the lesions become necrotic and crack. A whiteish ooze may exude from the cracks. The sample should be scraped locally from the rind of the fruit.
- **Bacterial samples:** Use a toothpick to remove a colony of bacteria from a culture plate. Stir into a microcentrifuge tube containing 300 µL of SEB4 extraction buffer. Do not use Aac ImmunoStrips with cell culture broth. Skip to step 5.

**Note:** If reusing cutting tools, disinfect them with a 10 % bleach solution between every sample.

**2.** Cut open the sample extraction bag near the bottom of the label. Be careful not to spill the buffer.  
 \*SEB4 buffer is required to perform this assay. (Figure 3)

**3.** Insert the sample between the mesh linings near the bottom of the sample extraction bag.

**4.** Extract the sample by thoroughly macerating it with an Agdia tissue homogenizer or a blunt object such as a pen or marker. (Figure 4)

An adequately extracted sample will result in a homogenous green or light brown colored solution.

**Allow the resulting solution to settle for 3 minutes before inserting the ImmunoStrip.**



Figure 1

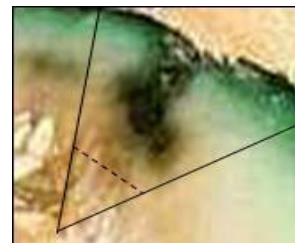


Figure 2



Figure 3



Figure 4

## Perform Assay

**5.** Remove an ImmunoStrip then reclose the container. When handling the ImmunoStrips, always grasp the top of the ImmunoStrip marked with Agdia's name. Do not remove the protective covering.

Insert sample end of the ImmunoStrip into the channel portion of the bag (no mesh) until submerged in the extract **up to the white line approximately ¼ inch from the bottom. Do not allow the side of the ImmunoStrip to come into contact with foam or bubbles (if present).** (Figure 4)

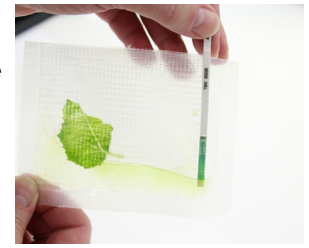


Figure 4

**6.** Place the bag in a letter holder or another device in an **upright** position. Allow the ImmunoStrip test to remain in the sample extract **for 30 minutes**. Positive results may be visible in as few as 5 minutes. Lower titer samples may take up to 30 minutes.

## Interpret Results

**7.** Remove the ImmunoStrip from the extract and interpret the results. Use the images provided as a guide to determine results. If storing the ImmunoStrips as a permanent record, immediately cut off the sample pad, then press the remaining ImmunoStrip between paper towels to remove any excess liquid.

If only the control line is visible, this indicates a negative result.

If the control line is visible and the test line is also present at any intensity of pink\*\*/purple, this indicates the presence of the target pathogen (or in some cases, a closely related pathogen). Visit the product webpage to see if any other pathogens are known to cross-react with this test.

The control line assures that the test is working properly. If the control line does not appear, the test is invalid, even if a test line is visible (see troubleshooting).

As with all diagnostic tools, Agdia recommends confirming all results with a secondary detection method before making any economic decisions (ex: discarding plants due to positive test results, etc.).

## SAFETY

Agdia recommends reading all relevant SDS sheets before using assay components: <http://docs.agdia.com/datasheets.aspx>.

## TROUBLESHOOTING

|  |   |
|--|---|
| Control line did not develop.            | <ul style="list-style-type: none"> <li>Submerging the ImmunoStrip past the white line in the sample extract. (Step 5)</li> <li>ImmunoStrip inserted before the 3 minute sample extract settling period. (Step 4)</li> </ul>   |
| Test runs very slow or not at all.       | <ul style="list-style-type: none"> <li>Extracting more tissue than is required. (Step 1) Further dilute sample extract 1:10 in SEB1 and repeat test.</li> <li>Components were not warmed to room temperature before use. (Kit Storage)</li> <li>Check kit and components expiration dates.</li> </ul>   |
| Test has a green or pigmented test line. | <ul style="list-style-type: none"> <li>Extracting more tissue than is required. (Step 1)</li> <li>Green lines should be considered a negative result. (Step 7)</li> <li>**In rare cases, red, orange, or purple fruits and tissues (for example, red cucumbers) may cause what appears to be a positive test line. Contact Agdia before testing these types of samples. (Step 7)</li> </ul> |
| Test and/or control line is weak.        | <ul style="list-style-type: none"> <li>Components absorbed moisture. (Kit Storage) Moisture can cause the membrane to wick without test components and fail to produce lines.</li> <li>Low pathogen titer in the sample. (Step 7)</li> <li>Check kit and components expiration dates.</li> </ul>  |

## QUESTIONS OR TECHNICAL SUPPORT:

Phone: 800-622-4342 (toll-free) or 574-264-2014

Fax: 574-264-2153

E-mail: [info@agdia.com](mailto:info@agdia.com) for sales and general product information  
[techsupport@agdia.com](mailto:techsupport@agdia.com) for technical information and troubleshooting

Web: [www.agdia.com](http://www.agdia.com)

ImmunoStrip® is a registered trademark of Agdia, Inc.

m225.2 Revised: 11/08/2022

Scan for How To Video



Page 2 of 2

## Positive Results    Negative Results    Invalid Results

